


Whiteness Tester C-130

Simple Guide to Sampling

This guide explains the steps to use for accurate sampling when using the C-130 Whiteness Tester. **Please be sure to read Section 6 "Measurement" on page 9 of the user manual before use.**

1 Collecting the sample

Heap the sample on the sample tray.

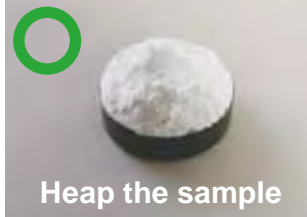

2 Packing down the sample.

Firmly press down on the sample using the included spatula.



If powder sticks to the outside or bottom of the sample tray, clean it with the brush.



Powder stuck to the sample tray may cause it to tilt, which will affect the reading.

3 Preparing the measurement surface.

In parallel motion, draw the spatula end toward you while firmly pressing down.



Contact with the air may cause drying or an increase in moisture content, thus changing the sample. As soon as sampling is complete, it is necessary to move quickly to the next phase of the operation (see page 10 of the user manual).